

DATA SHEET
Aqueous Mounting Medium
Cat. No. HK099-5K

Doc. No. HK099, Rev. No. D
Date of Release: 06-October, 2011

Reagents Supplied

HK099-5K	One bottle (15 ml) of Aqueous Mounting Medium, containing gelatin, glycerol and Sodium azide.
----------	---

Storage Conditions

Store at room temperature. For longer shelf life, store refrigerated.

Intended Use

This product may be used for aqueous mounting of frozen tissue sections, immunofluorescent preparations, or immunostains which require aqueous mounting. This product is a General Purpose Reagent suitable for diagnostic histopathology, laboratory and research use.

Method Of Use

1. Warm mounting medium for 10-15 minutes before use or until contents become liquefied. Do not exceed 75°C.
2. Remove slides from the last rinse with distilled water, and, without further drying, apply 2-4 drops of mounting medium to the tissue or cell specimen. Cover immediately with a coverslip.

General Purpose Reagent suitable for diagnostic histopathology, laboratory and research use.